
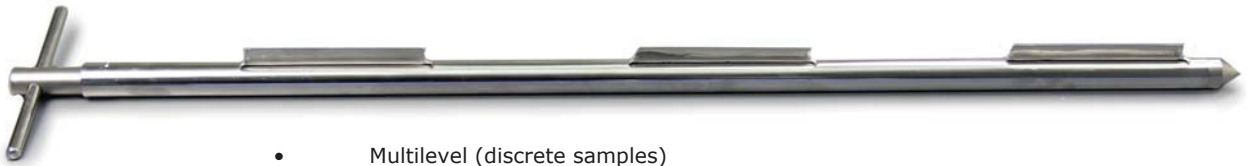


Cohesive Pocket Sampler

Zone sampling of Cohesive powders

The Cohesive Pocket Sampler has been specially adapted from the Pocket Sampler (on pages 36 & 37) to enable the taking of samples from damp and cohesive powders.

Scrapers have been added to push the sticky powder into the sample pockets. Scrapers are fully welded to give a hygienic, GMP correct sampler.

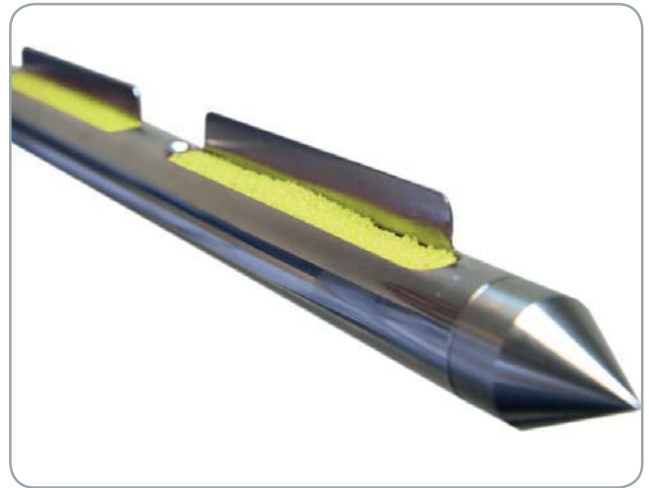


- Multilevel (discrete samples)
- 316 Stainless steel construction
- Suitable for sticky, cohesive powders

B900

Operation

1. Insert the sampler into the product
2. At the required depth rotate handle to expose the collecting pockets.
3. Rotate the sampler through 360° so the product is pushed into the collecting pockets
4. Rotate handle to close pockets and remove sampler.

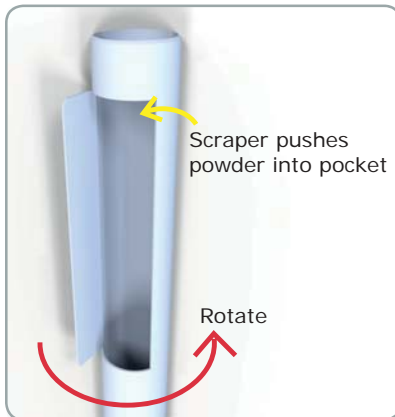

QUICK QUOTE

ONLINE PRICING

Stock No.	Length	No.Pockets	Pocket Volume	Material	Diameter
675 2220B600	600mm	3	15ml	316ss	25mm
675 2220B900	900mm	3	20ml	316ss	25mm
675 2220B1500	1500mm	5	20ml	316ss	25mm

675 TIP2 Rounded Tip (Sampler supplied with pointed tip)

Cleaning Brushes - see page 129

Sampler Cases - see page 130

XRF Certification - see page 139

* Diameter does not include scrapers. Scrapers are approximately 10mm high