



QAQC LAB

www.qclabequipment.com

SEED RIFFLE TYPE DIVIDER

The QAQC Lab RIFFLE DIVIDER is suitable for dividing dried seeds and grains into 2 equal halves, using two series of chutes.



Includes main body, two pans, and one scoop. Made from high quality stainless steel.

The scoop lays inside the hopper, and when removed, drops the grains on to the chutes.

Hopper Dimensions approximately 430 mm x 120 mm

Weight 15 lbs Packaged 15" x 22" x 18"

673 6718

Number of chutes 24

Width of chutes 12 mm

Hopper Size 430 mm x 120 mm

No of pans 2

Instructions: Place the scoop in the hopper in such a way that it is at a right angle from the hopper, by holding the handle of the scoop. The scoop will rest in the middle of the hopper.

Pour the sample into the hopper, and lift the scoop. The grain is divided into two equal divisions, and discharged into separate pans.



[QUICK QUOTE](#)

[ONLINE PRICING](#)